

PVT Next Day Legionella PCR™ Test Instructions

Next Day Legionella PCR™

IMPORTANT: Perform and ship this test **ONLY** on Monday, Tuesday, Wednesday, or Thursday

1 GATHER MATERIALS
One water sample location requires the following items found in the shipping carton:



250 mL Sterile Sample Bottle with Location ID Label



Security Seal



Shipping Label
(Fedex Priority Overnight)

In addition to the supplied items (above) you will need:


- Packing Tape Gloves
 Safety Glasses
(Recommended)

2 AFFIX LOCATION ID LABEL
Affix completed location ID label horizontally to water sample bottle.

3 COLLECT WATER SAMPLE
IMPORTANT: *Gloves are required and safety glasses are recommended to be worn.*
Fill water sample bottle to neck of bottle with water from source to be tested. Close bottle tightly and briefly shake (5-10 sec.) to dissolve Sodium Thiosulfate Neutralizing Powder.

4 APPLY SECURITY SEAL
Apply security seal to water sample bottle over side of cap to side of container exactly as shown. Make sure security seal does not cover location ID label.

5 REPACK and SHIP on SAME DAY
For assurance of safety in shipping, repack water sample bottle upright (lid up) in shipping carton as shown in the photo. Close and seal lid with tape. Apply Priority Overnight shipping label.

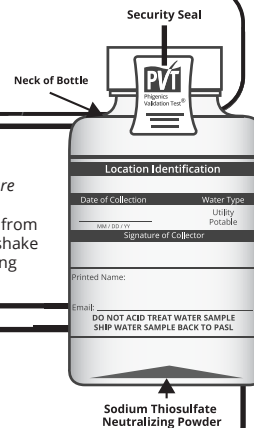

Neck of Bottle

Security Seal

Location Identification

Date of Collection Water Type
Utility
Portable

Printed Name:

Email: DO NOT ACID TREAT WATER SAMPLE
SHIP WATER SAMPLE BACK TO PAST

Sodium Thiosulfate
Neutralizing Powder

250 mL Water Sample Bottles