



Phigenics  
Validation Test®



## PVT Swab Collection Procedure

---

1. Remove the swab from the sterile tube.
2. Collect sample by rubbing the swab over the sample area, reversing directions between strokes. Repeat four more times over the same surface area.
3. Immediately replace the swab into the neutralizing buffer, making sure cap is tightly closed.
4. Affix tube with a completed location label.

### **5. Ship sample overnight to:**

Phigenics Analytical Services Laboratory  
35701 West Ave, Suite 100  
Warrenville, IL 60555