



# Copy your Reference Control Library to another computer for offline unmixing and analysis\*

\*Requires backup and restoration of Instrument Data/User Accounts to link associated QC values


**1** Login to **SpectroFlo** and go to **QC & Setup** module

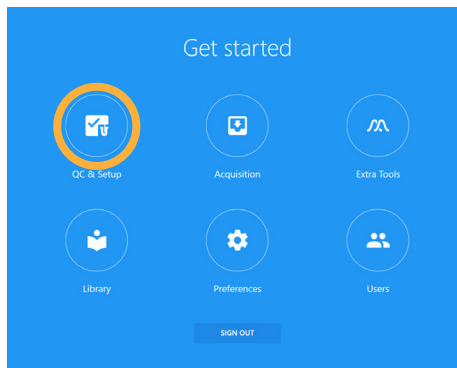


**2** Select **Reference Controls** then **Export**

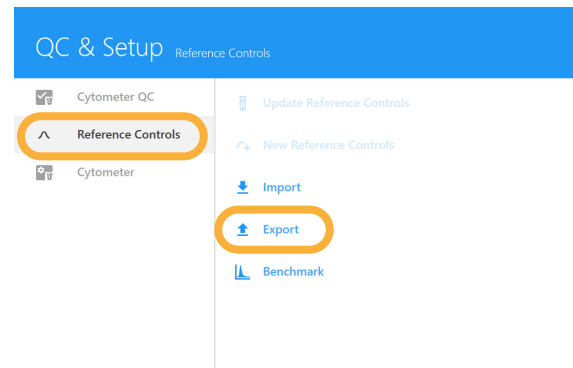


**3** Choose any (or all) controls you wish to export, **Save** the resultant zip file

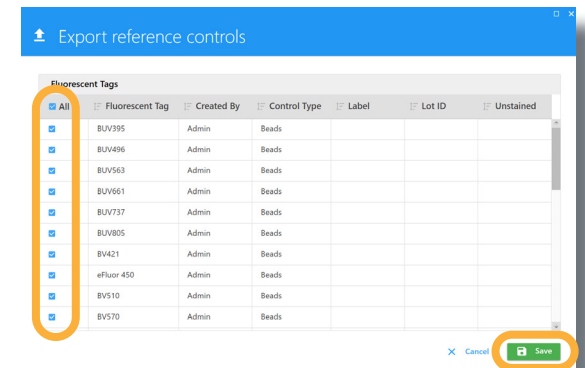

**4** At offline computer, login to **SpectroFlo** and go to **QC & Setup** module

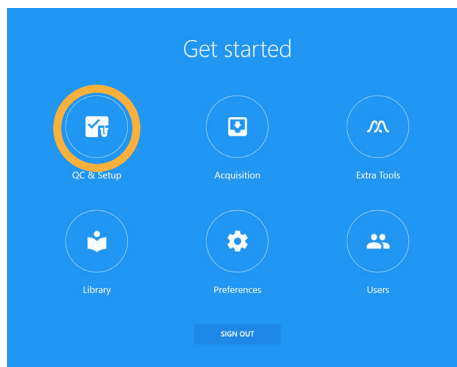

**5** Select **Reference Controls** then **Import**, select the zip file saved in step 3

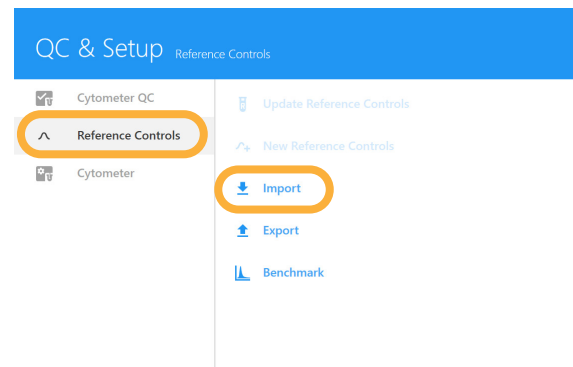

Get More Instructables  
cytekbio.com

