


# Praxis Moisture Balance

{ The User Guide

\*This guide is intended for use in addition to the Operating Instructions included with your Praxis Moisture Balance. It is not meant to replace the manufacturer's instructions, nor does it provide comprehensive instructions for the use of the Praxis Moisture Balance, aka the "Citizen Moisture Analyzer".

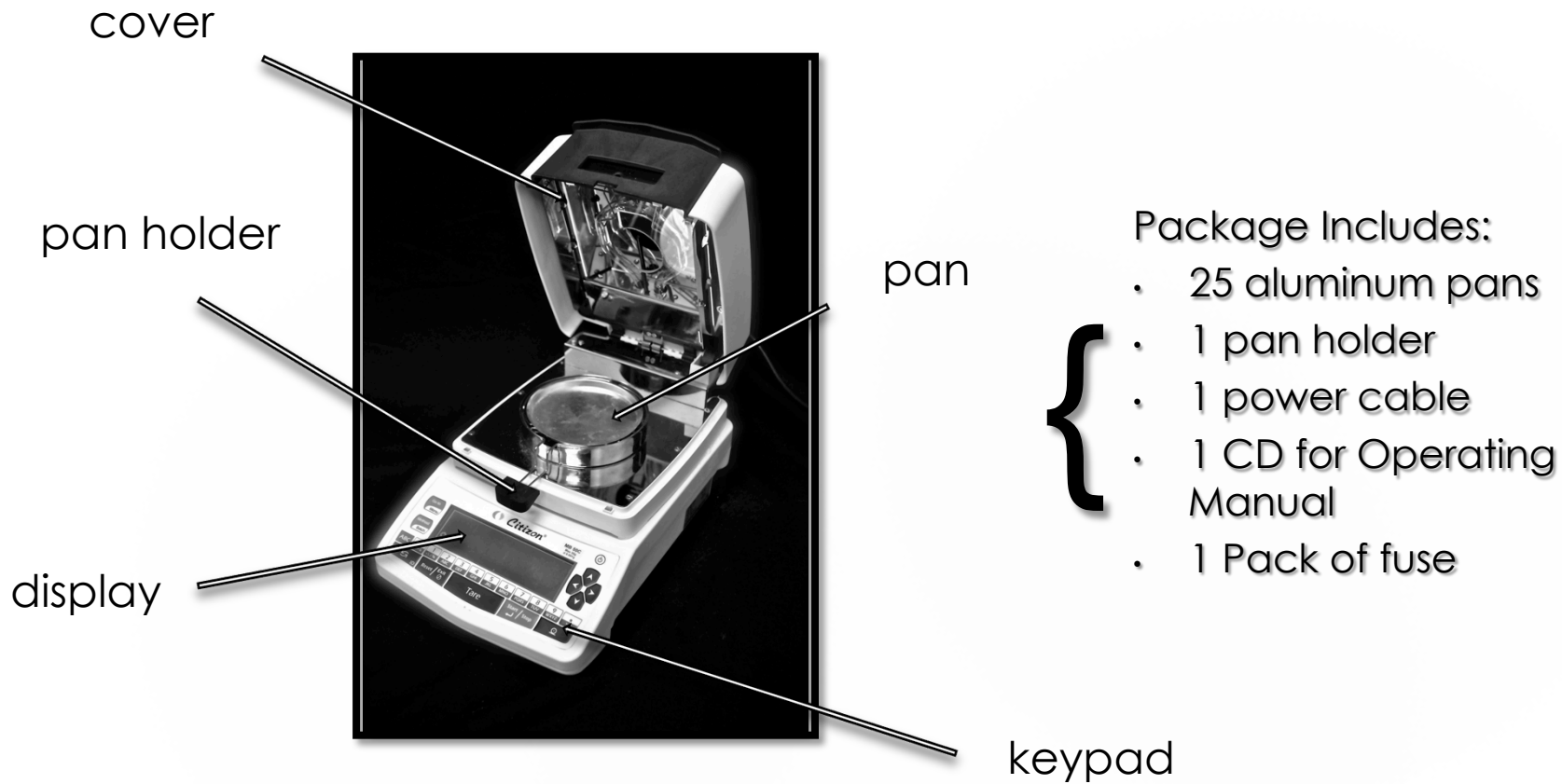

# The Basic Components

# The Buttons

{ Explained



- Wakes up analyzer
- Puts analyzer in stand-by mode

On/Off Key



- Press once for “Go-to” mode
  - Set functionality in Settings => Soft key
- Press and hold for two seconds for “Menu”
  - Displays four sub-menu choices:
    - Method
    - Settings
    - Data
    - Calibrate

# Go-to/Menu Key



- Press once for “Method” function
  - Displays method list
- Press and hold for two seconds for “Batch” function
  - Displays batch list
  - **Note:** data storage mode must be set to “Batchwise” to use this function

# Method/ Batch Key



- Activates letter/symbol mode or number mode
  - Keypad (highlighted in yellow) displays the text that the button will create in each mode
    - Black background is the letter/symbol mode
    - White background is the number mode

# ABC Key



- **Shown:** Keypad is in letter/symbol mode
- Pressing the ABC Key at this point would switch the keypad to number mode

Light background with dark letters indicates the active setting

# ABC Key






- In Pre-Analysis mode, press once to tare the weight of the pan
- Also used as the Backspace or Delete Key when typing with the keypad

# Tare Key



- Press to Start/Stop a test
- Also used as the Enter Key when typing with the keypad
- Press to select choices when changing the settings for your tests

# Start/Stop Key

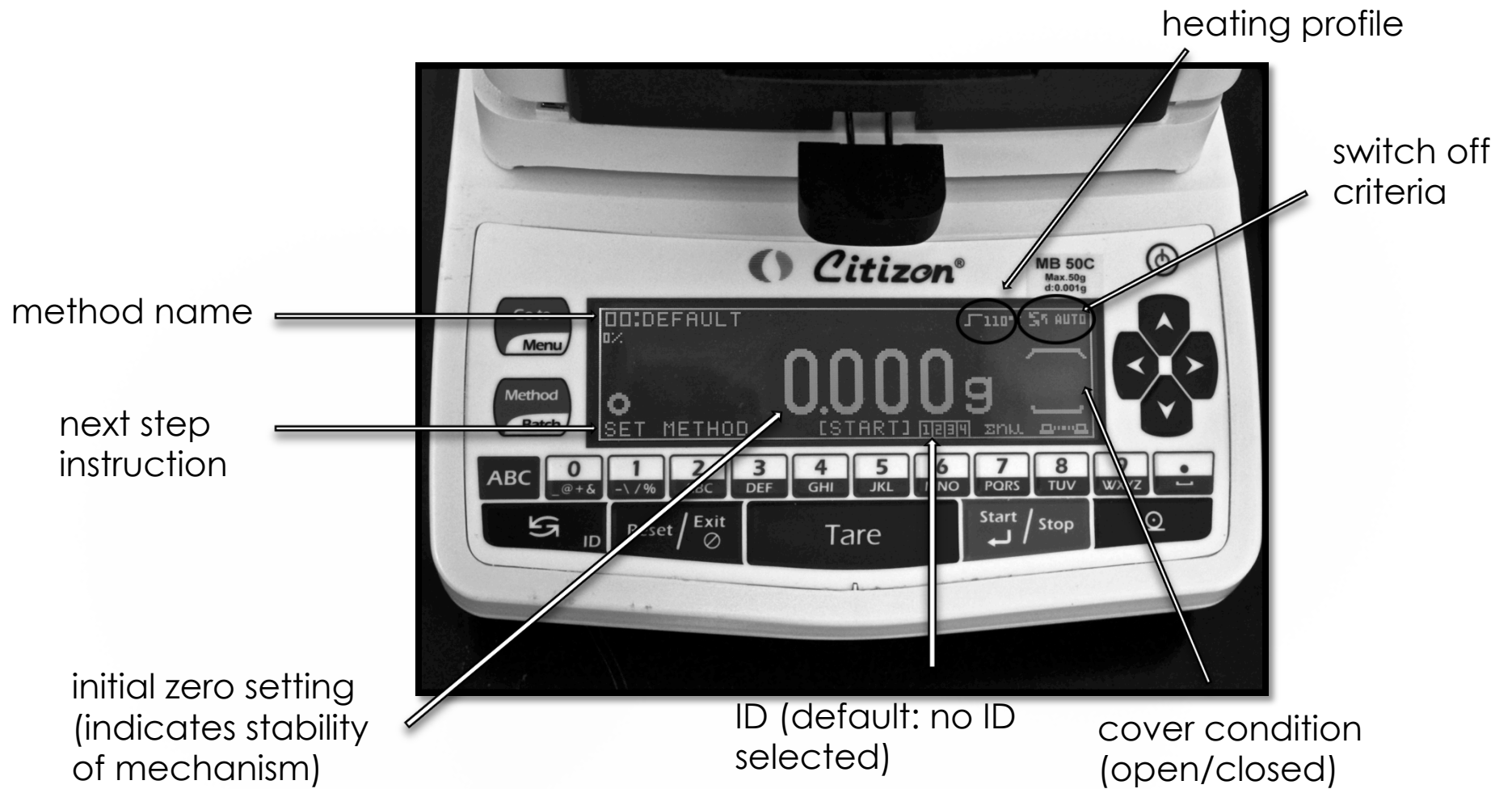

Shown: closed

# Initial Screen Display




- Press once to print results (must be connected to a printer via USB)

# Print Key

# Running a Test

{ Step-by-Step

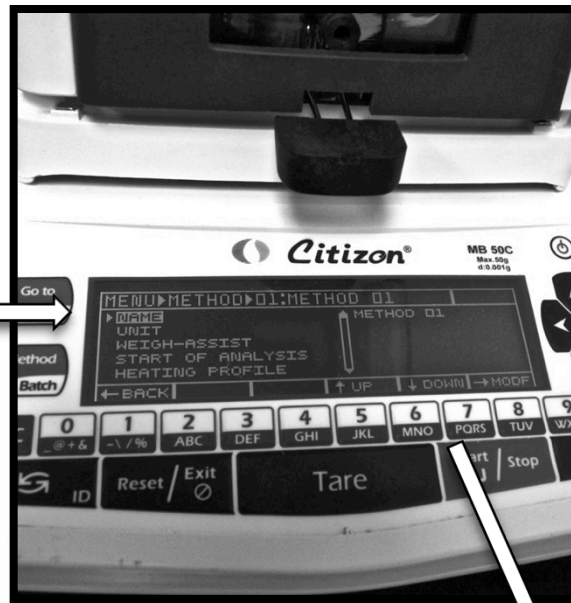

- Press "Menu"
- Choose "Methods"
- Select "Method 01" (or whichever number you wish to name)
- Select "Name"
- Delete the default method name by pressing "Tare" until the space is clear. Then, type in your choice of name using the keypad.
- Hit "Start/Stop" to save the name, then press the left arrow button to go back to set any other variables

# Step 1: Set the Method Name



### weigh-assist

- Set Target Weight
- Set Tolerance



### heating profile


- Standard
- Gentle
- Rapid
- Steps (multiple temperature settings during testing)



### standby heating

- On/Off

# Set Other Variables

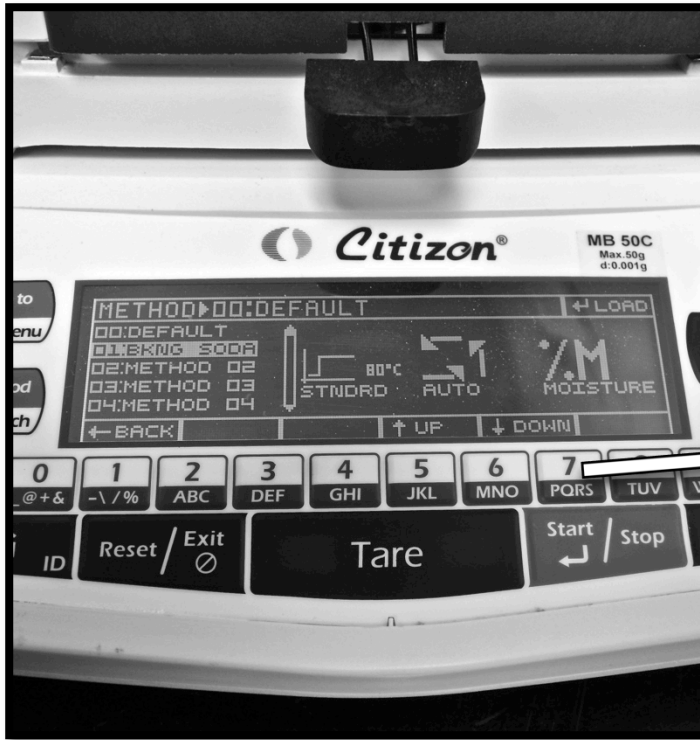

Press "Method". Then, use the arrow keys to move to your selection



Press "Start/Stop" to choose your selection. The new method will appear in the top left corner of the display screen

## Step 2: Choose the Method





Make sure to tare the empty pan weight before loading sample



Distribute the sample evenly on the pan

## Step 3: Prepare the Sample



{ Close the cover and press "Start/Stop".

## Step 4: Start the Test



Testing Complete!



{ Use a USB cable to connect the Praxis to a printer

## Step 5: Printing (part 1)



Press the "Print" button (circled) to print results

## Step 5: Printing (part 2)



Correct



Incorrect



After running the sample test, open the cover using the top lip (shown on left). Do not open using the pan holder.

**\*Helpful Hint\***